EN	ES	IT	NL
DA	NO	AR	JP
FR	DE	PT	EL
FI	SV	RU	UA
CS	HU	LV	PL
BG	SL	KOR	TH
ET	LT	MT	RO
SK	TR	MS	ZH

Single Use Garment Type PB[6]

INSTRUCTIONS FOR USE

BioClean-D™ BioClean-C™

Single Use Garment Type PB[6]

This instruction for use note is to be used in combination with the specific information that is mentioned on each packaging enclosure.

USE OF GARMENT

BioClean partial body protection garments (Type PB[6]) complement the BioClean range of full body protection (Type 5/6) garments. These are single use garments designed to protect against limited liquid splashes and fine sprays. Type PB[6] garments have not been subjected to the whole suit test. They are also designed to provide product protection. Check the garment carefully to ensure there are no defects. Ensure the expiry date has not been reached. For full protection wear compatible coverall, hood, gloves, facemask, goggles and footwear as appropriate.

WEARING GARMENT

- **Hood:** Pull fully over head. Ensure elasticated facial opening creates a good seal. For full protection wear compatible coverall, gloves, facemask, goggles and footwear offering at least equivalent protection. These should be taped together using waterproof tape to cover each join.
- **Overboot:** Place over foot and pull up leg. Where overboot has ties, secure to ensure comfortable fit. For full protection wear compatible hood, coverall, gloves, facemask and goggles offering at least equivalent protection. These should be taped together using waterproof tape to cover each join.
- **Sleeve cover:** Don through cuffed end where garment label is attached. Pull onto lower arm and ensure there is a good seal at wrist and above elbow. For full protection wear compatible hood, coverall, gloves, facemask, goggles and footwear offering at least equivalent protection. These should be taped together using waterproof tape to cover each join.
- **Apron:** Don apron (if applicable put arms into sleeves) and then secure at back of the neck using Velcro neck fastening, using tie-tapes at the waist tie securely at the rear.
- **Lab Coat:** Don garment and fasten studs.

LIMITATIONS OF USE:

Not suitable for prolonged or intensive exposure to chemical spray and splash. Heat stress should not occur, but consideration should be given to alternation of work and rest periods. The user is the sole judge of the suitability of the garment, and any accessories, given its intended use and duration. Nitritex Ltd accepts no responsibility for improper use.

GARMENT FABRIC:

Polyethylene/Polypropylene laminate.

WARNING (BioClean-D only):

The person wearing the electrostatic dissipative protective clothing shall be properly earthed. The resistance between the person and the earth shall be less than 10⁵ Ω.

Electrostatic dissipative protective clothing shall not be open or removed whilst in presence of flammable or explosive atmospheres or while handling flammable or explosive substances;

Electrostatic dissipative protective clothing shall not be used in oxygen enriched atmospheres without prior approval of the responsible safety engineer; The electrostatic dissipative performance of the electrostatic dissipative protective clothing can be affected by wear and tear, laundering and possible contamination.

Electrostatic dissipative protective clothing shall permanently cover all non-complying materials during normal use (including bending and movements).

STORAGE & TRANSPORT:

Store in a clean, dry place away from heat and light. To prevent the risk of damage, transport in cardboard box provided and retain unopened until required.

DISPOSAL:

Remove carefully and turn inside out to trap contamination. Assess any contamination that may have been introduced during use and destroy immediately in accordance with regulations.


To the best knowledge of the manufacturer, BioClean-D or BioClean-C material does not contain any constituent substances which could lead to skin irritations or skin diseases or other allergic reactions or that are carcinogenic, toxic to reproduction or mutagenous.

CAUTION:

Keep plastic bag away from children to prevent the danger of suffocation.

EXPLANATION OF GARMENT LABEL - KEY:

1. Coverall Manufacturer
2. Trademark
3. Model Identification
4. CE marking - complies with Regulation (EU) 2016/425 relating to Category III PPE. Under the supervision of Notified body - SGS Fimko Oy, P.O. Box 30 (Särkiniementie 3), 00211 HELSINKI, Finland (EU Notified Body No. 0598).
5. Wearer should read these instructions
6. Indicates compliance with EU standards for protective clothing: Type PB[6] standard EN 13034:2005+A1:2009 for reduced spray, limited splash & EN1149-5:2008 antistatic (BioClean-D only)
7. International Care Symbols a) Do not wash b) Do not chlorine bleach c) Do not tumble dry d) Do not iron e) Do not dry clean f) Inflammable – keep away from naked flame and intensive heat
8. Garment size
9. DO NOT REUSE

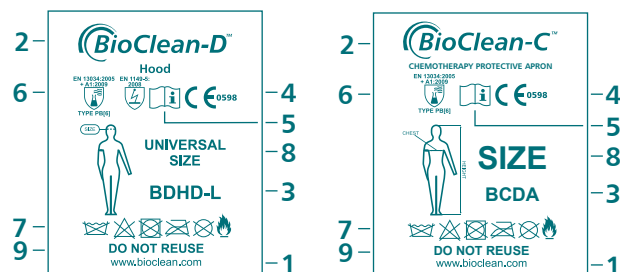The EU Declaration of conformity may be downloaded from:
www.ansell.com/regulatory

EU Type examination carried out by:
Centexbel-Gent, Technologiepark 70, BE-9052 Zwijnaarde (Gent), Belgium.
Notified Body Number 0493

TESTING

The **BioClean-D** range of PPE Category III Type 5/6 & Type PB[6] Single use garments have been subject to the following comprehensive tests:

Fabric Property Tests	Test Method	Class or Result
Abrasion Resistance	EN 530 Method 2	1
Trapezoidal Tear Resistance machine dir./x dir.	ISO 9073-4	3/2 (55.5N/29.3N)
Flex Cracking Resistance	ISO 7854 Method B	3
Seam Strength	ISO 13935-2	2 (70N)
Puncture Resistance	EN 863	1 (8N)
Tensile Strength Machine dir./x dir.	EN ISO 13934-1:1999	2/1 (97N/48N)
Charge Decay Time	EN 1149-3: 2004	0.07 sec
Particle Shedding	TEST-RP-CC-003.4 2011 (Helmke Drum)	1

Resistance to Penetration by Liquids EN 6530:2005

Liquid	Repellency Index	Penetration Index
Sulphuric Acid (30%)	3	3
Sodium Hydroxide (10%)	3	3
o-Xylene	2	3
Butan-1-ol	3	3

The **BioClean-C** range of PPE Category III Type PB[6] Single use garments have been subject to the following comprehensive tests:

Fabric Property Tests	Test Method	Class or Result
Abrasion Resistance	EN 530 Method 2	1
Trapezoidal Tear Resistance machine dir./x dir.	ISO 9073-4	3/2 (50.6N/32.3N)
Seam Strength	ISO 13935-2	2 (74N)
Puncture Resistance	EN 863	1 (10N)
Tensile Strength Machine dir./x dir.	EN ISO 13934-1:1999	2/1 (94N/47N)
Particle Shedding	TEST-RP-CC-003.4 2011 (Helmke Drum)	1
pH	ISO 3071: 2005	9.7
Azo Dyes	EN 14362-1: 2017	Pass

Resistance to Penetration by Liquids EN 6530:2005

Liquid	Repellency Index	Penetration Index
Sulphuric Acid (30%)	3	3
Sodium Hydroxide (10%)	3	3
o-Xylene	1	3
Butan-1-ol	3	3

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Ansell Healthcare Products LLC
Ansell Commercial México S.A. de C.V.